

Agilent D1K ScreenTape System Quick Guide

Principles

The Agilent 2200 TapeStation system is a tape-based platform for simpler, faster and more reliable electrophoresis. It is made up of three elements:

- 2200 TapeStation System (p/n G2964AA) or 2200 TapeStation Nucleic Acid System (p/n G2965AA)
- High Sensitivity D1K ScreenTape ~(p/n~5067-5363) with High Sensitivity D1K Reagents (Ladder and sample buffer) (p/n~5067-5364) or

D1K ScreenTape $(p/n \ 5067-5361)$ with D1K Reagents (Ladder and sample buffer) $(p/n \ 5067-5362)$

Agilent 2200 TapeStation Software

Kits

The High Sensitivity D1K ScreenTape (p/n 5067-5363) and D1K ScreenTape (p/n 5067-5361) are designed for analysing DNA from 35 – 1000 bp.

Specifications

Analytical Specification	High Sensitivity D1K ScreenTape	D1K ScreenTape	
Sizing Range	35 – 1000 bp	35 – 1000 bp	
Resolution ¹	35 – 300 bp: 15 %, 300 – 1000 bp: 10 %	35 – 300 bp: 15 %, 300 – 1000 bp: 10 %	
Sensitivity ²	5 pg/µL ³	0.05 ng/µL	
Sizing Precision	5 % CV	5 % CV	
Sizing Accuracy ⁴	±10 %	±10 %	
Quantitative Precision	15 % CV	10 % CV	
Quantitative Accuracy	±20 %	±20 %	
Quantitative Range	75 – 1000 pg/μL	0.1 – 50 ng/µL	
Carry Over	N/A	N/A	
Physical Specification			
Analysis Time	16 samples < 20 min, 96 samples < 100 min	16 samples < 20 min, 96 samples < 100 min	
Samples per consumable	16	16	
Sample Volume Required	2 μL	1 μL	
Shelf Life	4 months	4 months	
Box/Kit size	112 samples/box	112 samples/box	

 1 $\,$ Resolution is defined as the separation of fragments at half peak height or better

² signal:noise ratio > 3 for single peak

³ 2200 TapeStation Nucleic Acid System (G2965AA)

⁴ Determined using the D1K ladder as sample



Agilent D1K ScreenTape System Quick Guide

Essential Measurement Practices

Required tips and tubes for the TapeStation	 Optical Cap 8x Strip, Box of 120, 0.2 mL (p/n 401425) and Optical Tube 8x Strip, Box of 120, 0.2 mL (p/n 401428) Loading tips, 1 x384 (p/n 5067-5153) 		
Steps before use on the TapeStation	Equilibrate each vial to room temperature.Vortex mix each vial and briefly spin.		
Steps during sample preparation	• Keep reagents at room temperature during sample preparation.		
Storage after use on the TapeStation	 Store all reagent vials and ScreenTape at 2 – 8 °C Never store reagents and ScreenTape at room temperature or below 0 °C. If you run less than 16 lanes, store used tape upright at 2 – 8 °C for maximum of 2 weeks. 		
Pipette carefully	 Always pipette reagents against the side of the sample tube. If using a standard pipette ensure that no residual material is left on the outside of the tip. 		
Mix properly after each pipetting step	 Mix = Vortex the PCR tubes or 96 well plate on half-speed for 5 s. Spin = Move the samples to the bottom of the tubes/wells by pulsing in a centrifuge. 		

Storage Conditions

- Reagents vials: 2 8 °C
- ScreenTape: 2 8 °C (if you run less than 16 lanes, store used tape upright at 2 8 °C for a maximum of 2 weeks, never freeze ScreenTape any ScreenTape that is accidentally frozen should be discarded)

Products for analysing DNA

ScreenTape D1K and Reagents		
High Sensitivity ScreenTape D1K		7 ScreenTape
High Sensitivity D1K Reagents		1 vial
 High Sensitivity D1K Ladder 		75 μL
High Sensitivity D1K Sample Buffer		360 µL
and Reagents		
ScreenTape D1K		7 ScreenTape
D1K Reagents	1 vial	
D1K Ladder		75 μL
D1K Sample Buffer		360 µL
	High Sensitivity ScreenTape D1K High Sensitivity D1K Reagents • High Sensitivity D1K Ladder • High Sensitivity D1K Sample Buffer and Reagents ScreenTape D1K D1K Reagents • D1K Ladder	High Sensitivity ScreenTape D1K High Sensitivity D1K Reagents • High Sensitivity D1K Ladder • High Sensitivity D1K Sample Buffer and Reagents ScreenTape D1K D1K Reagents • D1K Ladder

Additional Consumables required for the 2200 TapeStation

- Loading tips, 10 x384 (p/n 5067-5152) / Loading tips, 1 x384 (p/n 5067-5153)
- Optical Tube 8x Strip, Box of 120, 0.2 mL (p/n 401428) and Optical Cap 8x Strip, Box of 120, 0.2 mL (p/n 401425) or 96 -well Sample Plates, Pack of 10 plates (p/n 5067-5150) and 96 -well Plate Foil Seal, Pack of 100 foils (p/n 5067-5154)

Additional Material Required (Not Supplied)

- Volumetric pipette
- Vortex mixer
- Micro-centrifuge

Safety Information

WARNING

Toxic agents

The handling of solvents, samples and reagents can hold health and safety risks.

- → When using/handling the ScreenTape and working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing).
- Always follow good laboratory practices and adhere to the guidelines established in your laboratory.
- → Refer to product material safety datasheets for further information.
- → The volume of substances should be reduced to the minimum required for the analysis.

CAUTION

Damage to the 2200 TapeStation


→ Use only the recommended tips, tubes and plates within the 2200 TapeStation instrument.

DNA Sample Preparation (High Sensitivity D1K)

Parts required	p/n	Description
	5067-5363	High Sensitivity D1K ScreenTape
	5067-5364	High Sensitivity D1K Reagents (Ladder and sample buffer)

1 Prepare Ladder

- **a** Aliquot a minimum of 3 µL High Sensitivity D1K Ladder (-) into the first tube/well.
- **2** Mix 2 μL sample buffer (**b**) with 2 μL DNA sample by vortex for 5 s.
- **3** Spin down to position the sample at the bottom of the tube.

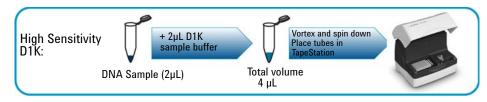


Agilent D1K ScreenTape System Quick Guide

DNA Sample Preparation (D1K)

Parts required	p/n	Description
	5067-5361	D1K ScreenTape
	5067-5362	D1K Reagents (Ladder and loading buffer)

- 1 Prepare Ladder
 - **a** Aliquot a minimum of 3 μL D1K Ladder (-) into the first tube/well.
- **2** Mix 3 μL sample buffer (**b**) with 1 μL DNA sample by vortex for 5 s.
- **3** Spin down to position the sample at the bottom of the tube.

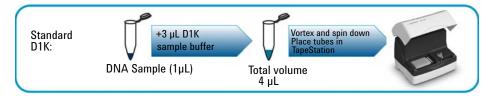

Sample Analysis

- 1 Launch the Agilent 2200 TapeStation software.
- 2 Load the samples, ScreenTape D1K and loading tips into the 2200 TapeStation.
- **3** Select the required samples on the controller software.
- 4 Click **Start** and specify a filename with which to save your results.

Limited Use Label License

Some products within this system contain SYBR® Green I, which is licensed from Molecular Probes Inc. for use in research and development only. SYBR® is a registered trademark of Molecular Probes, Inc.

Technical Support

For technical support, please visit www.agilent.com/genomics/contact.

Further Information

Visit Agilent Technologies` web site. It offers useful information, support and current developments about the products and technology: www.agilent.com/genomics/tapestation



Part Number: G2964-90030 Edition 10/11 **Printed in Germany** © Agilent Technologies, Inc. 2011 Agilent Technologies Hewlett-Packard-Straße 8 76337 Waldbronn, Germany