

Millennia ON/OFF Procedures (10W pumped Ti:S)

Su Lin's office #: 7-0391
Doug Daniel #: 5-4608

ON:

1. Turn on the chiller (located under the optical table, setting: 18C).
2. Press the 'Laser Power' button to warm up the diodes

The indicator will show the warming up progress

3. Hold the 'Laser Power' until you see the 'Laser Emission' on
4. Once you see this screen, use the up/Down button to set the power value (< 9.5W) and wait for the upper number to match the setting, reduce the power to 0.2 W if an error appears, and report the problem to me
5. Use the 'Setup' menu to open and close the shutter

OFF:

1. Close the shutter
2. Reduced the power back to 0.2 W
3. Press 'Laser Power' button to turn off the laser. The remote controller screen should be back to the screen shown on right
4. Turn off the Chiller

Millennia Remote Controller

